

Basic instructions to start run

1. `cd IN0DAQ` – Takes you to the DAQ program directory
2. `./InoC11` – Starts the DAQ run
3. After a little pause, you will be prompted for run number, please type run number, for ex. as 0026
4. Press any key to bring up the Menu
5. Press 3 (for enabling monitor interrupts) and return
6. Press 5 (for recording the data on disk) and return
7. Press 8 (for quitting from the menu) and return
8. You will get event data display on the upper part of the screen
9. After a gap of monitor run time (for ex. 100 secs), monitor data will be displayed on the lower part of the screen.
10. A sample screen shot with legends of various fields is given below.

Basic instructions to stop run

1. Press any key to bring up the menu
2. Press 7 to stop the run
3. Rename the temporary event data file to a regular file . For ex.
`mv temp.eve run0026.eve`
4. Rename the temporary monitor data file to a regular file. For
ex. `mv temp.mon run0026.mon`
5. You can use the following scripts to look at the event and
monitor data respectively.

```
OffDispEve <event data filename>  
OffDispMon <monitor data filename>
```